

How do I update my LabQuest original?

1. USB drive

- Download the current update software:
<https://www.scientrific.com.au/downloads.php?c=software>
- Get a fresh USB drive that is preferably no bigger than 2Gb (4Gb may be the smallest you can get now) and has **nothing else on it**.
- Place the LQ update files on the USB drive but ARRANGE them so that if your USB drive is drive D then the drive shows the files as:

D:\fw (a folder)

D:\fw_pre (a folder)

LQ_1.7.x_201x-yy-zz_Update.lqu (a file)

2. Battery

- Charge your LabQuest battery until it is fully charged before proceeding.
- If your battery is over 3 years old it might be time to purchase a new one.

3. Update procedure

- With the USB drive in the LabQuest's USB port turn the LQ on
- Tap on Home – Control Panel – System Information – Update LabQuest
- Select USB icon
- Scroll down to update file and select
- Ok – Yes to apply update
- Allow the update to proceed

4. Calibrate screen when prompted

5. Check update has been completed:

- Tap on Home – Control Panel – System Information
- Verify new update version number is displayed

6. Charge LabQuest overnight